

# NEBNext<sup>®</sup> for Ion Torrent<sup>™</sup>

LIBRARY PREPARATION KITS

# Table of Contents

3	General Introduction
4–5	DNA Library Preparation Workflow
6	Product Selection
6–7	Product Details
8–9	Additional NEBNext Features
10	Technical Tips
11	Ordering Information

## TOOLS & RESOURCES

Visit [NEBNext.com](https://www.nebnext.com) to find:

- The full list of products available
- Protocols & FAQs
- Online tutorials to help with product selection, general handling tips and more
- NEBNext citations



The NEBNext suite of products supports Ion Torrent sequencing with library preparation tools that streamline workflows, minimize inputs, and improve library yields and diversity. DNA library preparation is complete in approximately 2 hours, with minimal hands-on time. In addition to the extensive QCs on individual kit components, all NEBNext kits for Ion Torrent are functionally validated by preparation of a library, followed by Ion Torrent sequencing.

### NEBNext for Ion Torrent formats:

Sets include reagents for the entire library preparation workflow, and include master mix reagents that minimize the number of components and pipetting steps.

### Bulk & Custom formats:

Choose NEBNext's bulk and custom formats when your reagent needs exceed standard volumes, or you require a specialized formulation. As enzyme manufacturers, we are able to easily provide customized reagents, kits and modules to meet your specific NGS library prep workflow needs. For more information, please contact [NEBsolutions@neb.com](mailto:NEBsolutions@neb.com).


#### THE NEBNext ADVANTAGE

- Broad range of input amounts, from low ng to µg
- Fast workflows with minimal hands-on time
- Gel-free workflows
- High yields
- High library diversity
- Minimized GC bias
- Convenient formats include kits and modules
- All reagents undergo stringent quality controls, plus sequencing validation
- Value pricing



# Workflow for Ion Torrent DNA Library Preparation

Recommended input amounts	Total DNA
NEBNext Fast DNA Library Prep Set for Ion Torrent	10 ng – 1 µg
NEBNext Fast DNA Fragmentation & Library Prep Set for Ion Torrent	10 ng – 1 µg

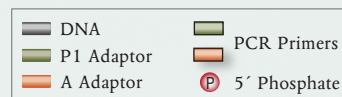


## 1 DNA Fragmentation

The first step in DNA library preparation is generally fragmentation of purified DNA. Fragmentation can be accomplished by a number of methods including acoustic shearing, nebulization and enzyme-based methods, such as dsDNA Fragmentase®. All of these methods leave a mix of 3' and 5' ends (recessed, overhang, blunt) which may or may not be phosphorylated. NEBNext kits for Ion Torrent are available with or without enzyme-based fragmentation reagents, allowing a choice of fragmentation method.




## 2 End Repair and 5' Phosphorylation

Blunt-ended fragments are created by filling in or chewing back 3' and 5' overhangs. Phosphorylation of the 5' ends ensures the fragments are suitable for ligation.

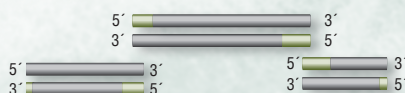

## 3 Adaptor Ligation and Nick Translation

Short adaptor sequences are added to the ends of the DNA fragments by blunt-end ligation. The adaptors are not 5' phosphorylated, in order to minimize adaptor-dimer formation. Since ligases require a 5' phosphate for ligation, this absence of 5' phosphate results in a nick on one strand at each ligation site. This nick is repaired by nick translation, using Bst 2.0 WarmStart™ DNA Polymerase. The four possible reaction products are shown below.

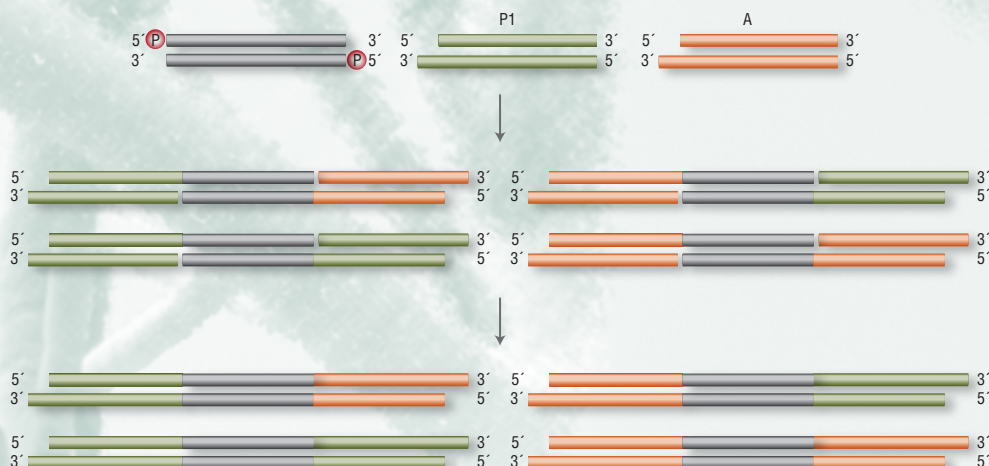


## 4 PCR Enrichment

Finally, amplification using a high-fidelity polymerase is performed, resulting in a robust library suitable for sequencing.

This step has multiple functions:

- Selection for molecules with the proper adaptor at each end
- Increase the amount of library
- Incorporation of sequences required downstream in the sequencing workflow, and, when desired, barcodes/indices to enable multiplexing. NEBNext kits for Ion Torrent are not supplied with barcode adaptors, but are compatible with adaptors from other sources, including Ion Xpress™ Barcode Adaptors from Life Technologies.

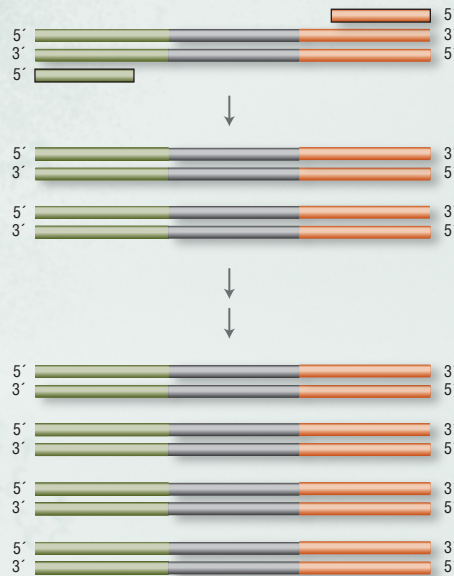

## Product Selection

Designed with the user in mind, NEBNext kits maximize efficiency and convenience. For use with Ion Torrent, NEBNext kits are available with or without fragmentation reagents, allowing the user to choose the method of fragmentation. For maximum convenience and ease of scale-up, the NEBNext Fast DNA Fragmentation & Library Prep Set for Ion Torrent (NEB #E6285) includes a master mix for enzyme-based fragmentation. This set contains dsDNA Fragmentase, a mix of two enzymes for random DNA shearing, described on page 9, and this is combined with End Repair reagents for a streamlined workflow. When mechanical shearing methods, such as Covaris® acoustic shearing or nebulization are preferred, the NEBNext Fast DNA Library Prep Set for Ion Torrent (NEB #E6270) is the recommended choice. The components and protocol for this kit have been optimized for the volumes and concentrations of DNA sheared by mechanical methods. Both kits have the same fast workflow and low minimum input DNA requirements.

## Product Details

In addition to stringent QC on individual components, the NEBNext DNA kits are functionally validated by library preparation of a genomic DNA library, followed by Ion Torrent sequencing. Reagent lots are reserved specifically for inclusion in NEBNext kits. Most of these reagents are provided in master mix format, reducing the number of vials provided in the kits, and reducing pipetting steps. Adaptors and primers for singleplex libraries are supplied in the kits. For multiplexed libraries, the Ion Xpress Barcode Adaptors from Life Technologies can be used.



**Input Amount**  
10 ng – 1 µg\*

Fragmentation/End Repair

Adaptor Ligation/Fill-In

PCR Enrichment

NEBNext Fast DNA Fragmentation & Library Prep Set for Ion Torrent (NEB #E6285)		
<b>Library Prep Set</b>	<ul style="list-style-type: none"> <li>DNA Fragmentation Master Mix</li> <li>DNA Fragmentation Reaction Buffer</li> </ul>	<ul style="list-style-type: none"> <li>Adaptors for Ion Torrent</li> <li>T4 DNA Ligase</li> <li>T4 DNA Ligase Buffer for Ion Torrent (10X)</li> <li><i>Bst</i> 2.0 WarmStart DNA Polymerase</li> </ul>
		<ul style="list-style-type: none"> <li>Primers for Ion Torrent</li> <li>NEBNext High-Fidelity 2X PCR Master Mix</li> </ul>



**Hands-On Time**  
12 min.  
**Total Time**  
110 min. – 133 min.

NEBNext® Fast DNA Library Prep Set for Ion Torrent™ (NEB #E6270)		
<b>Library Prep Set</b>	<ul style="list-style-type: none"> <li>End Repair Enzyme Mix</li> <li>End Repair Reaction Buffer (10X)</li> </ul> <p>Reagents for Fragmentation are not included</p>	<ul style="list-style-type: none"> <li>Adaptors for Ion Torrent</li> <li>T4 DNA Ligase</li> <li>T4 DNA Ligase Buffer for Ion Torrent (10X)</li> <li><i>Bst</i> 2.0 WarmStart DNA Polymerase</li> </ul>
		<ul style="list-style-type: none"> <li>Primers for Ion Torrent</li> <li>NEBNext High-Fidelity 2X PCR Master Mix</li> </ul>



**Hands-On Time**  
12 min.  
**Total Time**  
110 min. – 133 min.

\* Note that a minimum of 100 ng is recommended when used in conjunction with Ion Xpress Barcode Adaptors.

Library Preparation workflow for Ion Torrent is complete in approximately 2 hours



**Input Amount**



**Time**

Fragmentation/  
End Repair/  
Heat Inactivation

Adaptor  
Ligation/  
Nick Translation

Clean Up/  
Size Selection

Amplification  
(Optional)


Clean Up



**Workflow Time**


10 ng – 1 µg	<b>Hands-On</b>					<b>Hands-On</b>
	2 min.	1 min.	5 min.	1 min.	3 min.	12 min.
	<b>Total</b>					<b>Total</b>
	30 min.	20 min.	44 min.	0 – 23 min.	16 min.	110 – 133 min.

Sequencing Data

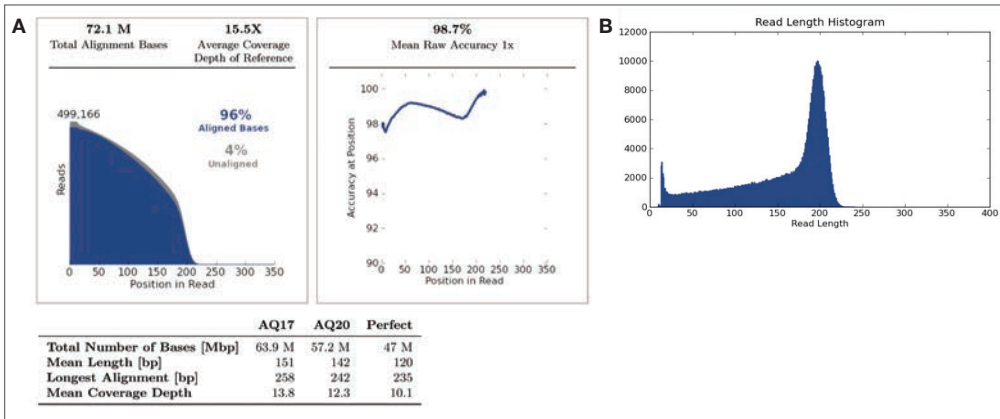

A typical Ion Torrent run report for libraries made from *E. coli* (K12 MG1655 strain) genomic DNA using NEBNext Fast DNA Fragmentation & Library Prep Set for Ion Torrent. A. Alignment Summary. 72 Mb of data was generated, with an average genome coverage of 15.5 X, from approximately 0.5 Million 200 bp reads. B. Read length histogram from 200 bp run.

Varying GC Content Libraries

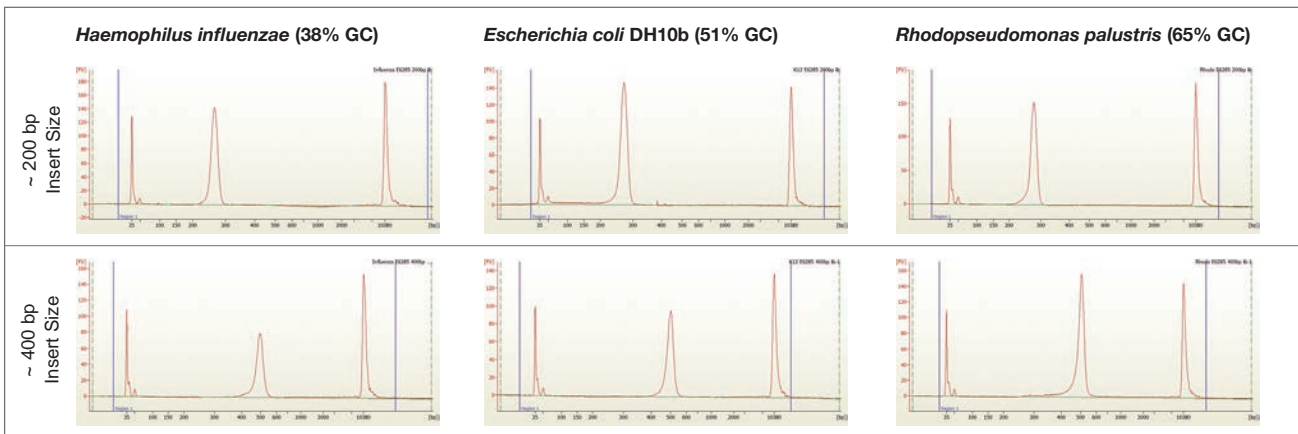


0.5 mg of DNA from 3 different genomes with varying GC content were used to construct 200 bp and 400 bp libraries using the NEBNext Fast DNA Fragmentation and Library Prep Set for Ion Torrent, analyzed by the Agilent® Bioanalyzer®.

## Minimized GC bias with the NEBNext Formulation of Q5<sup>®</sup> High-Fidelity DNA Polymerase


To ensure that sequence data reflects exactly the sequence of the original sample, it is essential that amplification of libraries be performed evenly and with high fidelity. Historically, high-fidelity polymerases have been more susceptible to difficulties in PCR amplification of GC-rich and other challenging regions. If such bias occurs in library amplification, this can lead to uneven sequence coverage, challenges in sequence assembly and even “missing” sequence.

To minimize GC bias, NEB offers the NEBNext High-Fidelity 2X PCR Master Mix (NEB #M0541). This master mix is specifically optimized for robust, high-fidelity amplification of next-generation sequencing libraries, regardless of GC content. This unique formulation combines the ultra low error rates of Q5 High-Fidelity DNA Polymerase with specific buffers and reaction enhancers to support robust amplification of a broad range of templates, even those with high GC content. The buffer component of the master mix has been optimized for robust, even amplification of NGS libraries, including GC-rich amplicons. This combination makes NEBNext High-Fidelity 2X PCR Master Mix an ideal choice for your NGS library construction.

### NEBNext High-Fidelity 2X PCR Master Mix minimizes GC bias and error rate

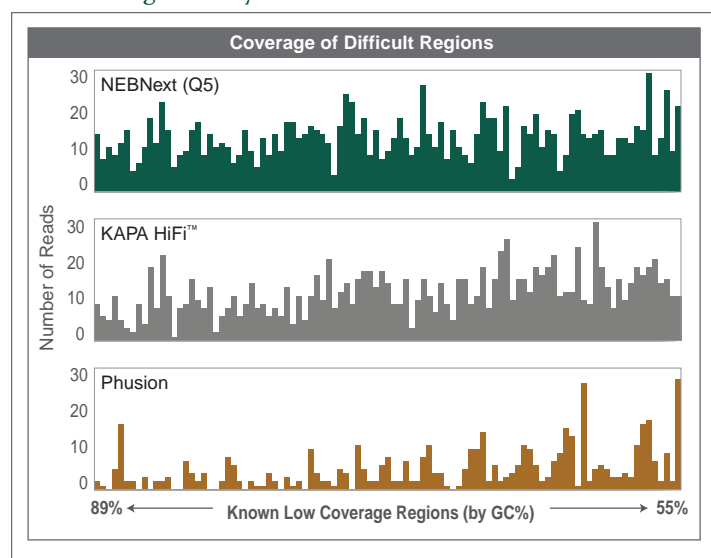

Indexed human IMR90 DNA libraries were split and amplified using 8 cycles of PCR with Phusion High-Fidelity PCR Master Mix with HF™ Buffer, KAPA HiFi<sup>®</sup> HotStart ReadyMix or NEBNext High-Fidelity 2X PCR Master Mix, followed by sequencing on an Illumina<sup>®</sup> HiSeq™ 2000. 180 million reads were randomly extracted from each data set and the number of reads overlapping distinct low coverage regions of the human genome (Aird et.al. *Genome Biology*, 2011) are shown for each library.

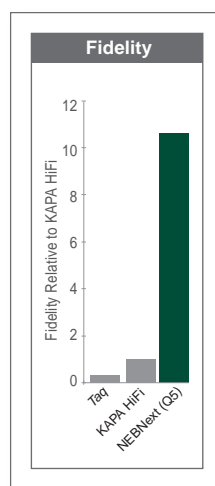

Fidelity measurements of Taq DNA Polymerase (in Standard Taq Buffer), KAPA HiFi HotStart ReadyMix and NEBNext High-Fidelity 2X PCR Master Mix were measured, side-by-side, in a PCR-based mutation screening assay using a lacZ method. Values ( $n \geq 2$ ) are expressed relative to KAPA HiFi HotStart ReadyMix.




## DNA Fragmentation with NEBNext dsDNA Fragmentase

Methods for DNA fragmentation include mechanical shearing (such as acoustic shearing or nebulization) and enzyme-based fragmentation.


For this step in the DNA library preparation workflow, NEB provides NEBNext dsDNA Fragmentase, an enzyme-based reagent that generates fragments of the desired sizes for NGS libraries in a time-dependent manner. dsDNA Fragmentase contains two enzymes; one randomly generates nicks on dsDNA, and the other recognizes the nicked sites and cuts the opposite DNA strand across from the nick, producing dsDNA breaks. The resulting DNA fragments contain short overhangs, 5'-phosphate and 3'-hydroxyl groups. Randomness of shearing for this application is critical, and dsDNA Fragmentase provides random fragmentation, similar to mechanical methods (1,2).

### Fragmentation of HeLa cell gDNA using dsDNA Fragmentase

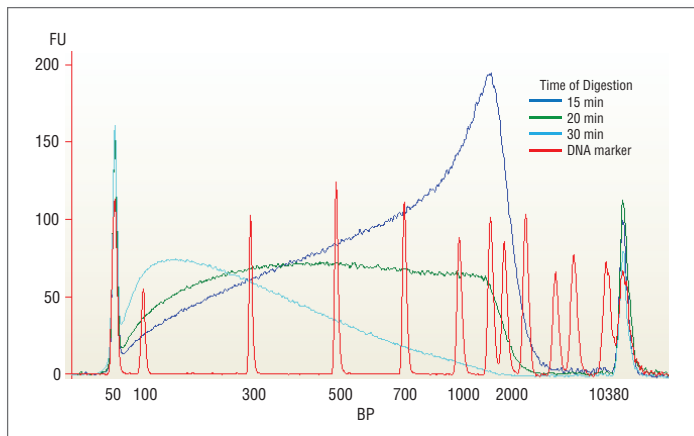

5  $\mu$ g of genomic HeLa DNA was incubated with NEBNext dsDNA Fragmentase for varying times as indicated. Fragmented DNA was purified using MinElute<sup>®</sup> columns and analyzed by the Agilent Bioanalyzer 2100.

### NEBNext dsDNA Fragmentase generates fragments in the 100–300 bp range more effectively than nebulization

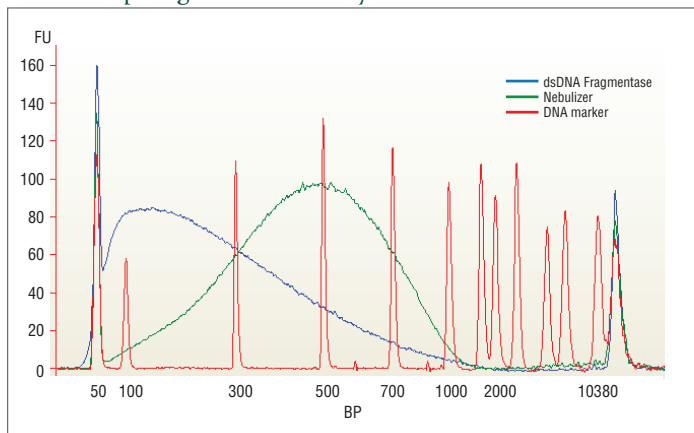

Relative size distribution of *E. coli* DNA fragments with dsDNA Fragmentase vs. nebulization as seen using the Agilent Bioanalyzer. The dsDNA Fragmentase sample was incubated for 30 minutes at 37°C with 0.05  $\mu$ g of DNA per  $\mu$ l of dsDNA Fragmentase in 1X Fragmentase Reaction Buffer with 100  $\mu$ g/ml BSA. The Nebulizer sample was prepared by nebulization of DNA in 50% glycerol for 6 minutes at 35 psi.

#### References

1. Kuirim, E. et al. (2011) PLOS One 6, e28240.
2. Adey, A. et al. (2010) Genome Biology 11: R119.

## General Handling Tips

- Spin down the vials before opening. This ensures that everything is at the bottom of the vial and not in the lid or on the sides.
- Use filter tips
- Use RNase- and DNase-free plastics, water, and solutions
- Thaw and keep all master mixes on ice
- Keep all sample and reagent vials capped when not in use
- Mix reaction components by pipetting gently up and down, and do not vortex (except when using Fragmentase, which should be vortexed as described in the protocol)
- Return enzymes to the freezer after use
- Always use freshly prepared 80% ethanol
- Keep all sample and reagent vials capped when not in use
- Always use aseptic technique.
- Store final libraries in TE, in low-bind DNA tubes

## DNA Sample Input Guidelines

### Integrity of DNA

- Start with as high quality DNA as possible. The quality of the input material directly affects the quality of the library. DNA quality can be assessed using gel electrophoresis. Impurities such as detergents or probes, or damaged DNA can usually be seen as a smear on the gel. RNA impurities are often seen at the bottom of the gel. Absorbance measurements can also be used as an indication of DNA purity. Ideally, the ratio of the absorbance at 260 nm to 280 nm should be between 1.8 – 2.0. However, measurements can be affected by the presence of RNA or small nucleic acid fragments.

### Quantitation of DNA

- It is important to accurately quantify the DNA sample prior to library construction. Fluorescence based detection which utilizes dsDNA specific dyes, such as the Qubit® from Life Technologies, are more accurate than UV spectrometer based measurements, as the presence of RNA or other contaminants can result in overestimation of the DNA sample.

### Bead-based clean-ups and size selection

- Be careful when transferring material not to disturb the bead pellet
- Be sure to vortex the beads just before use – they should be a uniform suspension
- Do not over-dry the beads. This can make resuspension difficult and reduce yield.

### TOOLS & RESOURCES

#### Visit NEBNext.com to:

- View the latest videos from NEB scientists discussing technical tips for library preparation



## Ordering Information

KITS FOR ION TORRENT DNA LIBRARY PREPARATION		NEB #	SIZE
DNA	NEBNext Fast DNA Library Prep Set for Ion Torrent	E6270S/L	10/50 rxns
	NEBNext Fast DNA Fragmentation & Library Prep Set for Ion Torrent	E6285S/L	10/50 rxns
MODULES & ENZYMES		NEB #	SIZE
DNA	NEBNext High-Fidelity 2X PCR Master Mix	M0541S/L	50/250 rxns
	NEBNext dsDNA Fragmentase	M0348S/L	50/250 rxns

FRAGMENTASE®, NEW ENGLAND BIOLABS® and NEBNEXT® are registered trademarks of New England Biolabs, Inc.

HF™ and WARMSTART™ are trademarks of New England Biolabs, Inc.

QUBIT® is a registered trademark owned by Life Technologies, Inc.

ION TORRENT™ and ION XPRESS™ are trademarks owned by Life Technologies, Inc.

AGILENT® and BIOANALYZER® are registered trademarks of Agilent Technologies.

COVARIS® is a registered trademark of Covaris.

MINELUTE® is a registered trademark of the Qiagen Group.

ILLUMINA® and HISEQ® are registered trademarks of Illumina, Inc.

KAPA HIFI® is a registered trademark of Kapa Biosystems.

## USA

New England Biolabs, Inc.  
Telephone (978) 927-5054  
Toll Free (USA Orders) 1-800-632-5227  
Toll Free (USA Tech) 1-800-632-7799  
Fax (978) 921-1350  
e-mail: [info@neb.com](mailto:info@neb.com)  
[www.neb.com](http://www.neb.com)

## Canada

New England Biolabs, Ltd.  
Toll Free: 1-800-387-1095  
[info.ca@neb.com](mailto:info.ca@neb.com)

## China, People's Republic

New England Biolabs (Beijing), Ltd.  
Telephone: 010-82378265/82378266  
[info@neb-china.com](mailto:info@neb-china.com)

## France

New England Biolabs France  
Telephone : 0800 100 632  
[info.fr@neb.com](mailto:info.fr@neb.com)

## Germany & Austria

New England Biolabs GmbH  
Free Call: 0800/246 5227 (Germany)  
Free Call: 00800/246 52277 (Austria)  
[info.de@neb.com](mailto:info.de@neb.com)

## Japan

New England Biolabs Japan, Inc.  
Telephone: +81 (0)3 5669 6191  
[info@neb-japan.com](mailto:info@neb-japan.com)

## Singapore

New England Biolabs, PTE. Ltd.  
Telephone +65 6776 0903  
[sales.sg@neb.com](mailto:sales.sg@neb.com)

## United Kingdom

New England Biolabs (UK), Ltd.  
Call Free: 0800 318486  
[info.uk@neb.com](mailto:info.uk@neb.com)

[www.neb.com](http://www.neb.com)

